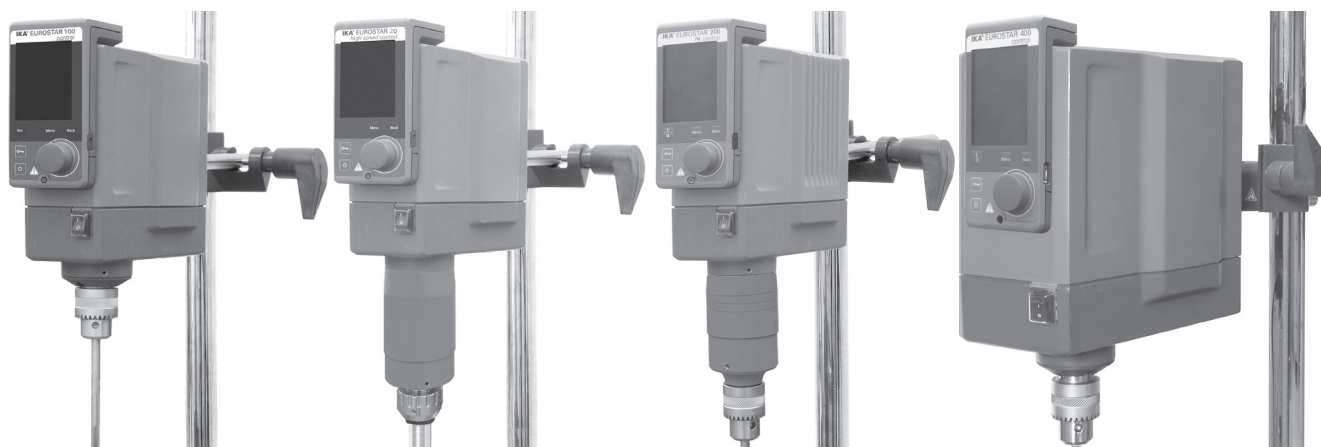
20000003848

EUROSTAR control_062015

IKA®

IKA® EUROSTAR 60 control
IKA® EUROSTAR 100 control
IKA® EUROSTAR 200 control
IKA® EUROSTAR 200 P4 control
IKA® EUROSTAR 20 high speed control
IKA® EUROSTAR 400 control

Kurzanleitung Ursprungssprache	DE	4			
User guide	EN	9	Kezelési útmutató	HU	64
Guide de l'utilisateur	FR	15	Navodila za uporabo	SL	68
Руководство пользователя	RU	19	Návod na obsluhu	SK	72
Guía del usuario	ES	24	Kasutusjuhend	ET	76
Istruzioni d'uso	IT	28	Lietotāja rokasgrāmata	LV	80
Gebruikershandleiding	NL	32	Vartotojo instrukcija	LT	84
Manual	SV	36	ръководство на потребителя	BG	88
Brugervejledning	DA	40	Manual de utilizare	RO	93
Bruksanvisning	NO	44	Οδηγίες χρήσης	EL	97
Käyttöohjeet	FI	48	用戶指南	ZH	102
Guia do usuário	PT	52	사용자 가이드	KO	106
Instrukcja	PL	56	ユーザーガイド	JA	111
Návod k obsluze	CS	60			

IKA®-Werke, Germany
Reg. No. 004343

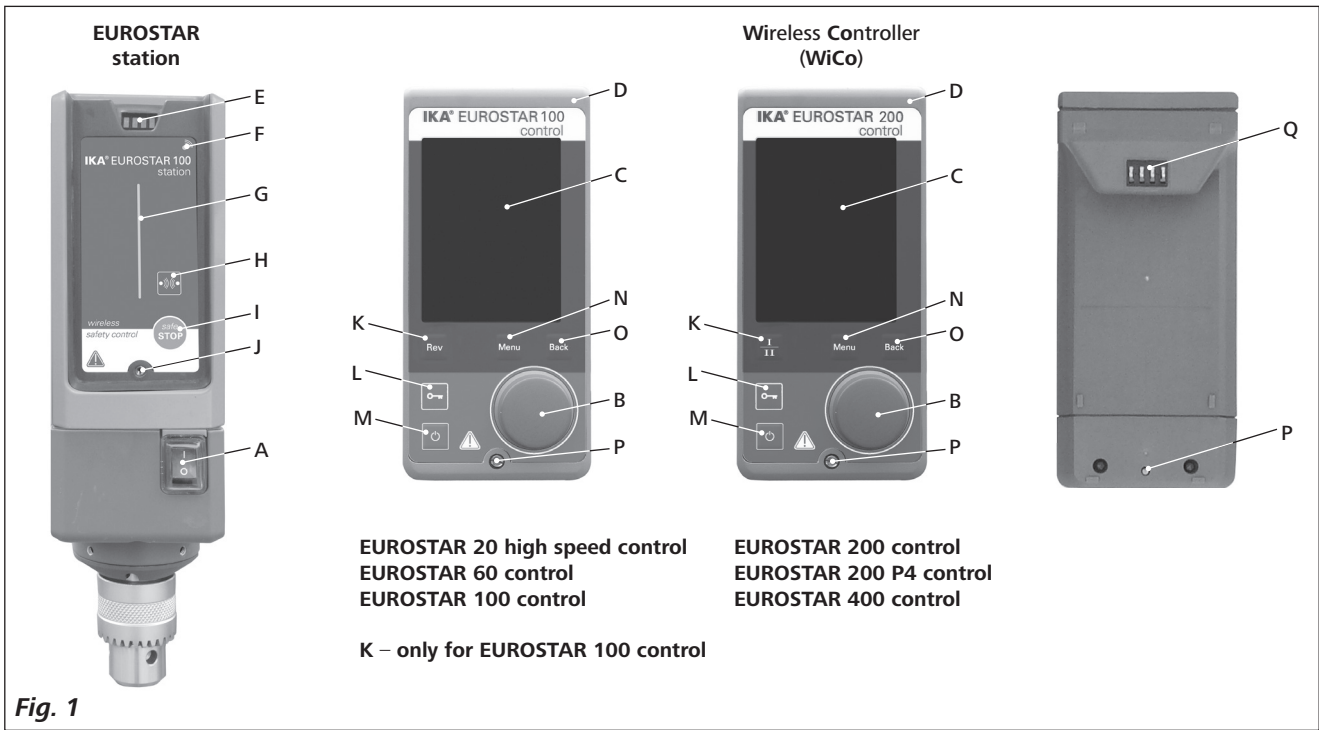
Fig. 1

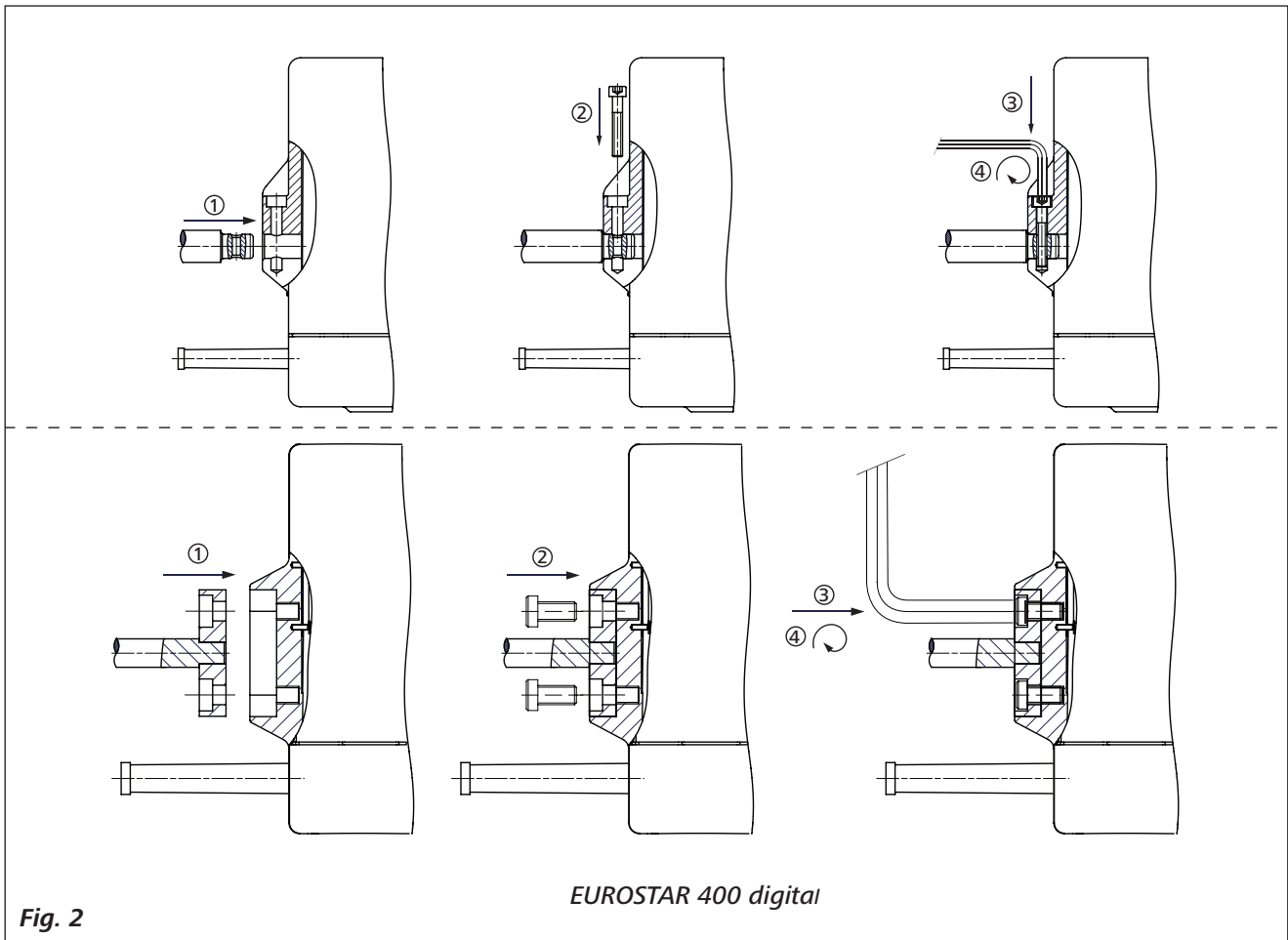
Fig. 2

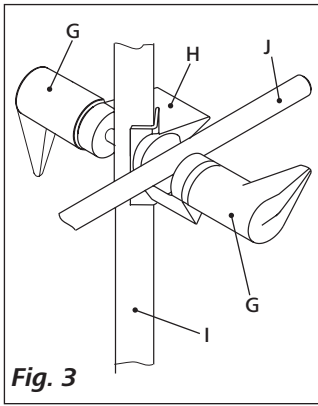
Fig. 3

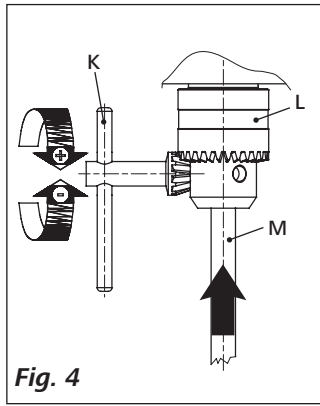
Fig. 4

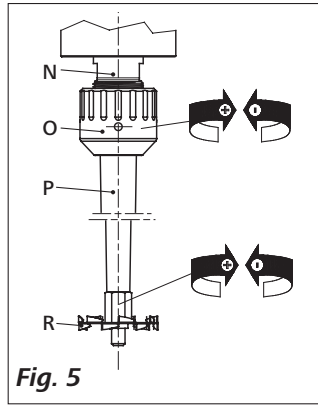
Fig. 5

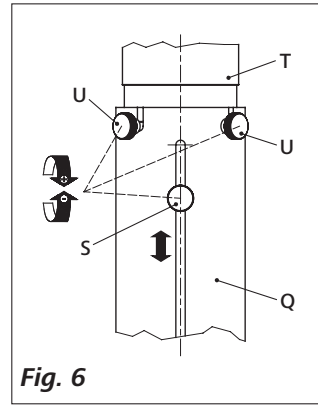
Fig. 6

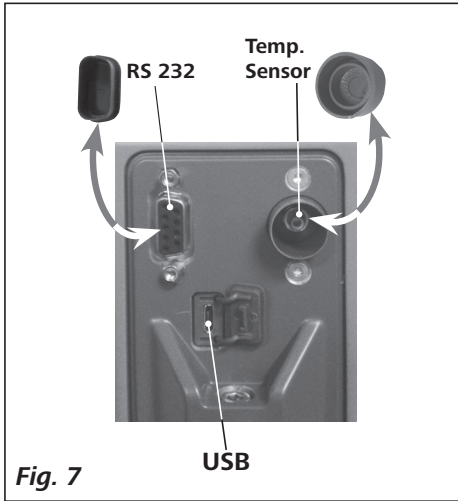
Fig. 7

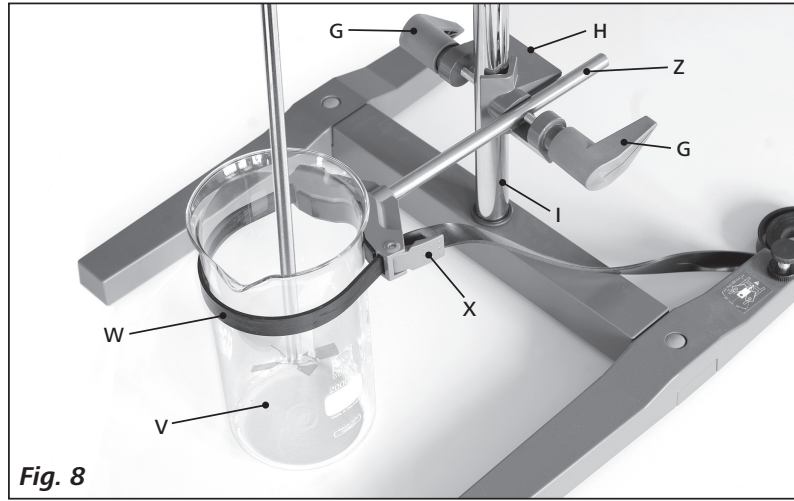
Fig. 8

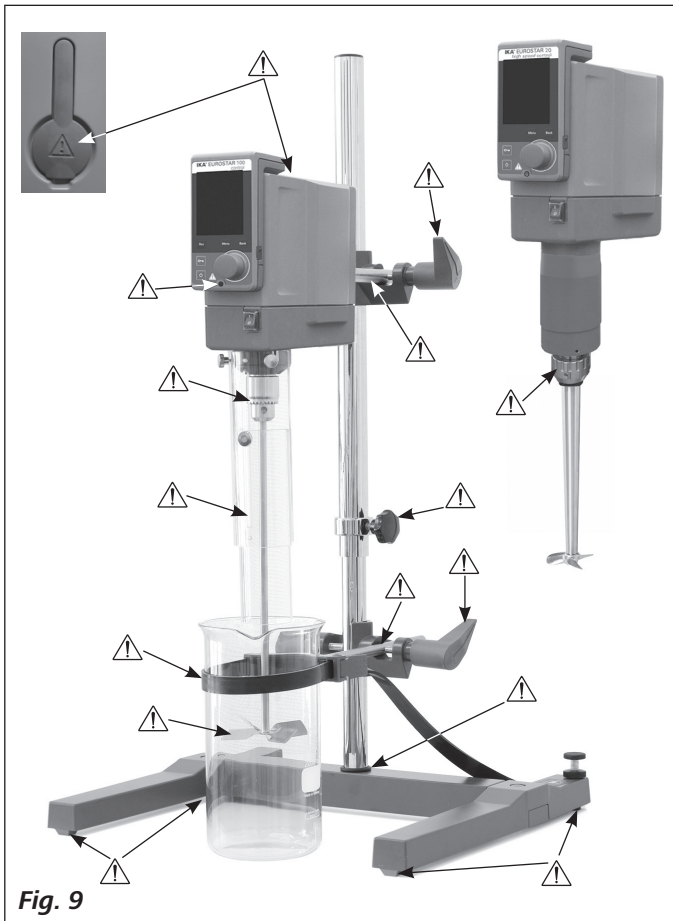
Fig. 9

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This user guide includes only the main operations and functions!

You can find and download **the complete operating instructions** in your language on our website (www.ika.com)!

Source language: German

	Page
EC Declaration of Conformity	9
Note for USA (FCC)	9
Note for Canada (IC)	9
Explication of warning symbols	10
Safety instructions	10
Correct use	12
Commissioning	12
Menu structure	14

EC Declaration of Conformity

We declare under our sole responsibility that this product corresponds to the regulations 2006/42/EC and 2004/108/EC and conforms with the standards or standardized documents EN 61010-1, -2-051; EN ISO 12100-1, -2; EN 60204-1 and EN 61326-1. Bluetooth® module:
Directive: 1999 / 5 / EC

Standards: EN 60950-1, EN 300 328, EN 301 489-1, -17

Note for USA (FCC)

This equipment complies with Part 15 of the FCC rules. Any changes or modifications not expressly approved by the Manufacturer could void the user's authority to operate the equipment. This device complies with Part 15 of the FCC rules subject to the following two conditions:

- (1) This device may not cause harmful interference, and
- (2) This device must accept all interference received, including interference that may cause undesired operation.

NOTE:

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment.

This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Note for Canada (IC)

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions:

- (1) This device may not cause interference
- (2) This device must accept any interference, including interference that may cause undesired operation of the device.

This device complies with Health Canada's Safety Code 6 / IC RSS-210. The installer of this device should ensure that RF radiation is not emitted in excess of the Health Canada's requirement.

Explication of warning symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the instrument or to system components.

This symbol indicates information **which is important for ensuring that the operations of the instrument are performed efficiently and for using the instrument**. Failure to observe this information can result in inaccurate results.

Safety instructions

- **Read the operating instructions in full before starting up and follow the safety instructions.**

- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the instrument.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Because the options for combining products, tools, stirring vessel, experiment and medium are nearly endless, user safety cannot be ensured simply with design requirements on the part of the product. For this reason, it may become necessary for users to take other precautionary safety measures. For example, glass device or other stirring vessels that are sensitive to mechanical stress can be damaged or shattered by an imbalance, increasing the speed too quickly or too little distance between the stirring element and the stirring vessel. Users can suffer serious injury from glass breakage or from the freely rotating stirring element.
- Uncontrolled reactions can be triggered by mixing the heated material insufficiently or by the energy generated by selecting a speed that is too high. In case of these and other increased operational hazards, users must take additional appropriate safety precautions (e.g. shatter protection). In any case, when using critical or hazardous materials in your processes, **IKA®** recommends to use additional appropriate measures to ensure safety in the experiment. For example, users can implement measures that inhibit fire or explosions or comprehensive monitoring equipment. Furthermore, users must make sure that the **OFF** switch of the **IKA®** product can be accessed immediately, directly and without risk at any time.

If installation or positioning cannot ensure this access at all times, an additional **EMERGENCY STOP** switch that can be easily accessed must be installed in the work area.

- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- Do not operate the instrument in explosive atmospheres, with hazardous substances or under water.
- Process pathogenic materials only in closed vessels under a suitable fume hood. Please contact **IKA®** application support if you have any question.
- The instrument is not suitable for manual operation.

- The high torque developed by the **EUROSTAR** requires particular care in the choice of stand, cross sleeve and anti-rotation element for the agitating vessel.
- Set up the stand in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- You must ensure that the stirring element is securely clamped in the chuck!
- Use stirring shaft protective equipment!
- The agitated vessels used for stirring have to be secured. Consider on a good stability of the entire structure.

Please pay attention to the dangerous parts of the equipment in **Fig. 9**.

- Protect the instrument and accessories from bumping and impacting.
- Check the instrument and accessories beforehand for damage each time when you use them. Do not use damaged components.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" section.
- Always switch the main switch in the **OFF** position or disconnect the power before changing stirring element and fitting allowed accessories.
- The instrument can only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
- The socket for the mains cord must be easily accessible.
- Socket must be earthed (protective ground contact).
- The voltage stated on the type plate must correspond to the mains voltage.
- Please observe the permitted speed for the stirring element. Never set higher speed.
- Make certain that the unit is set at the lowest speed before commissioning; otherwise, the unit will begin running at the speed set in last operation. Gradually increase the speed.
- Pay attention when setting the speed to any imbalance of the stirrer tools and possible spraying of the medium to be stirred.

Never operate the instrument with the stirrer tools rotating freely. Ensure that parts of the body, hair, jewelry or items of clothing cannot be trapped by the rotating parts.

The operation of a free rotating shaft end is dangerous. Therefore, for safety reason, only insert through the stirring tool over the upper edge of housing at standstill.

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed, otherwise there is a risk of:

- splashing of liquids
- projectile parts
- body parts, hair, clothing and jewelry getting caught.

Beware of the risk of:

- flammable materials
- glass breakage as a result of mechanical shaking power.

Reduce the speed if:

- the medium splashes out of the vessel because the speed is too high
- the instrument is not running smoothly
- the instrument begins to move around because of dynamic forces
- an error occurs.

Do not touch rotating parts during operation!

- There may be electrostatic activity between the medium and the output shaft which could cause a direct danger.
- After an interruption in the power supply or a mechanical interruption during a stirring process, the unit does not restart automatically.
- It is important to note that the surfaces of the motor (cooling fins) and certain parts of the bearing may get very hot during operation.
- Never cover the ventilation slots or cooling fins on the motor or on the instrument.
- Avoid knocking and impacting on the lower end of the shaft and the chuck gear teeth. Even minor, invisible damage can lead to imbalance and uneven shaft action.
- Ensure that the stand does not start to move.
- Imbalance of the output shaft, the chuck and in particular the stirring tools can lead to uncontrolled resonant vibrational behavior of the instrument and the whole assembly. Glass apparatus and stirrer containers can be damaged or shattered by this. It can cause injury to the operator, also can damage the rotating stirring tool. In this case exchange the stirring tool for one without imbalance or remedy the cause of the imbalance. If there is still imbalance, return it to the dealer or the manufacturer along with a description of the fault.
- If the instrument is operated too long in overload or if the ambient temperature is too high, the instrument switches off permanently.
- The machine must only be opened by trained specialists, even during repair. The instrument must be unplugged from the power supply before opening. Live parts inside the instrument may still be live for some time after unplugging from the power supply.

Covering or parts that are capable of being removed from the unit without accessory equipment have to be reattached to the unit for safe operation in order to prevent, for example, the ingress of fluids, foreign matter, etc..

If during operation the **battery pack RB 1** (rechargeable battery) becomes fully discharged, the device will continue to run or is shut down depending on the value settings for exceeding the time and safety speed. If the device is set so that it continues to run when the battery of the **Wireless Controller (WiCo)** is fully discharged, the only means of switching the station off are the "**safe STOP**" and the off switch!

Please note the following safety instructions for the battery pack RB 1 (rechargeable battery):

- Keep the battery pack out of reach of children at all times.
- Store the battery pack in a cool, dry place.
- Never throw the battery pack into a fire. Keep it away from direct sunlight and temperatures above 60 °C. High temperatures will damage the battery pack and render it unusable. Temperatures above 100 °C may cause it to explode.
- Never throw the battery pack into water or expose it to moisture. Water may lead to a short-circuit, causing the battery pack to explode.
- Do not deform or crush the battery pack or damage it in any other way. This can cause battery fluid to leak and/or the battery pack to explode.
- When not in use, keep battery packs away from paperclips, coins, keys, nails, screws or other small metal objects which could cause the contacts to be bridged. Short-circuiting may result in an explosion.
- Explosion of a battery pack may release battery fluid and cause a fire.
- The lithium polymer battery pack must only be used and charged in **IKA®** products designed for use with this battery pack.
- When the battery pack is inserted it should slide in easily and without resistance. Do not force it.
- If the battery pack is removed for an extended period of time, store it in a sealed plastic bag to prevent short-circuiting due to moisture or contact with metal components.
- The operating temperature range of the battery pack is from 0 °C to +45 °C. Note that the battery pack capacity will be reduced at temperatures below 20 °C.
- Only the rechargeable battery types recommended in the technical data may be used in the device!

Do not charge batteries that have leaked or that are discolored, deformed or damaged in any other way.

Disposal instructions:

- When disposing of the **IKA®** battery pack, please tape over the contacts with adhesive tape to prevent short-circuiting due to moisture or contact with metal components. Short-circuiting may result in an explosion.
- Do not throw used battery packs into your household waste. Dispose of them properly in accordance with statutory regulations.
- End users are obliged by law to return all used disposable and rechargeable batteries. Throwing them into the household waste is prohibited. Disposable/rechargeable batteries containing harmful substances are marked with this symbol to indicate that they may not be disposed of as household waste.
- You can return used disposable and rechargeable batteries free of charge to your local authority collection site or to any battery retailer. In doing so you will be complying with statutory regulations and helping to protect the environment.
- Batteries must be disposed of in accordance with local and national regulations.

Correct use

- **Use:**

For mixing/stirring liquids with low and high viscosities by various stirring tools.

Intended use: Stand device (chuck points down).

- **Range of use (indoor use only):**

- Laboratories
- Schools
- Pharmacies
- Universities

- **Wireless remote control:**

Before using the wireless link between the **Wireless Controller (WiCo)** and the laboratory device, first check whether your region is included in the radio communications approval for the device. If it is not, remote control can also be performed using a USB cable.

This instrument is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the instrument is operated with accessories that are not supplied or recommended by the manufacturer
- If the instrument is operated improperly or contrary to the manufacturer's specifications
- if the instrument or the printed circuit board are modified by third parties.

Commissioning

Assemble the overhead stirrer and all necessary accessories on a stable, even, non-slip surface. The **EUROSTAR** overhead stirrer must be secured to a stable stand (e.g. **R 2722** or **R 2723**) with a cross-sleeve (e.g. **R 270**). The stirring vessel must always be securely fixed for safety reasons. You must also ensure that the mounting device (stand) is set up in such a way that it is not liable to topple and does not start to move during the stirring procedure. Accessories must be assembled according to the assembly instructions (**Fig. 2** to **Fig. 8**).

Check whether the voltage given on the type plate corresponds to the available mains voltage.

The socket used must be earthed (fitted with earth contact).

If these conditions have been met, the machine is ready for operation when the mains plug is plugged in.

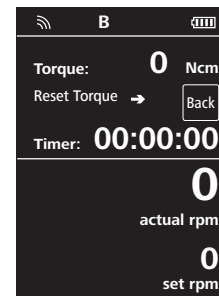
If these conditions are not met, safe operation is not guaranteed and the machine could be damaged.

Before the first time to use the **EUROSTAR** stirrer, fasten the **Wireless Controller (WiCo)** to the station with the screw (**P**, see **Fig. 1**), so that the rechargeable battery (**RB1 battery pack**) in the **Wireless Controller** can be charged.

After the device has been switched on at the main switch (**A**, see **Fig. 1**) the device name and software version appear in the display (**C**, see **Fig. 1**) of the **Wireless Controller (D**, see **Fig. 1**).

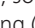
After This the following working screen appears automatically in the display.

Rotate the rotary/push knob (**B**, see **Fig. 1**) to set the target speed. Press the rotary/push knob (**B**), the stirrer starts to run with the set speed.

If the **Wireless Controller** is switched on and is not attached to the stirrer (the station), the green LED bar (**G**, see **Fig. 1**) and the green Bluetooth® LED (**F**, see **Fig. 1**) on the stirrer (station) will light up. Make sure that the speed selected is appropriate for the trial run. If in doubt, use the rotary knob (**B**, see **Fig. 1**) to select the slowest speed. To start or stop the stirrer function, press the rotary knob (**B**, see **Fig. 1**).

The controls of the **Wireless Controller** can be disabled by pressing the button (**L**) , so that no inadvertent changes are made while the device is running (the key symbol  appears in the display). Pressing the button (**L**)  a second time enables the controls again (key symbol  disappears from the display).

WARNING

In emergency the stirrer (station) can be switched off by pressing the "safe STOP" button (**I**, see **Fig. 1**) on the front of the stirrer. In this case the LED bar (**G**, see **Fig. 1**) changes color from green to red, and flashes.

A message appears in the display, advising that the **EUROSTAR** stirrer (station) has been forcibly switched off. To restart the stirrer, switch the main switch (**A**, see **Fig. 1**) of the **EUROSTAR** stirrer (station) OFF and ON.

If the Bluetooth® function of the **Wireless Controller** is active, the user can use the Bluetooth® search button (**H**)  to search for the **Wireless Controller**. A beep signal will be audible even if the **Wireless Controller** is switched off.

EUROSTAR 100 control:

The direction of rotation of the **EUROSTAR 100 control** stirrer can be changed by pressing the “Rev” button (**K**) when the stirrer is in stand-by status (stationary).

EUROSTAR 200, 200 P4 and 400 control:**• Speed setting:**

You can use the rotary knob (**B**, see **Fig. 1**) to pre-set the required speed before you start the device. If you then press the rotary knob (**B**, see **Fig. 1**) the device starts at the speed you set. When a change is made to the speed, the rated speed is shown in the display (**C**, see **Fig. 1**). When the device is stopped, the button (**K**, see **Fig. 1**) can be used to switch between the two speed ranges (**I** and **II**). When in standby mode, the display (**C**, see **Fig. 1**) shows the set speed.

The stirrer has two different speed ranges:

Range I: low speed, high torque.

Range II: high speed, low torque.

• Correct procedure for changing the speed range:

- Switch the device off at the rotary knob (**B**, see **Fig. 1**).
- Switch the speed range with button (**K**, see **Fig. 1**).
- Change the speed with the rotary knob (**B**, see **Fig. 1**).
- Switch the device on at the rotary knob (**B**, see **Fig. 1**).
- The speed can be varied at any time in operation.
- The speed appears in the display (**C**, see **Fig. 1**).

Menu structure

		Factory settings				
	Stirring	Speed Limit	Manual	2000 rpm EUROSTAR 60 control		
			Anchor stirrer	1300 rpm EUROSTAR 100 control		
			Centrifugal stirrer	2000 rpm EUROSTAR 200 control		
		Torque Limit	60 Ncm EUROSTAR 60 control			
			100 Ncm EUROSTAR 100 control			
		Torque Display	200 Ncm EUROSTAR 200 control			
			660 Ncm EUROSTAR 200 P4 control			
		Intermittent Mode	Run/Stop	20 Ncm EUROSTAR 20 high speed control		
			CW/CCW ❶	400 Ncm EUROSTAR 400 control		
			Interval	activated		
		Torque Calibration	Run Time	00:00 [mm:ss]		
			Stop Time	00:00 [mm:ss]		
			Temperature	Probe Temperature	-	
				Display	-	
			Timer	Set	00:00:00 [hh:mm:ss]	
				Display	activated	
			Operating Mode	A	-	
				B	activated	
				C	-	
			Display	Torque	activated	
Temperature	-					
Timer	activated					
	Programs	Program 1	-			
		...	-			
		Program 5	-			
	Safety	Time Out	00:30 [mm:ss]			
		Safe Speed	50 rpm EUROSTAR 200 P4 control			
		Password	100 rpm the other EUROSTAR			
		Vibration Sensor ❷	000			
	Settings	Languages	English	activated		
			Deutsch	-		
		Units	Temperature	°C	activated	
			°F	-		
		Display	Background ground	Black	activated	
				White	-	
			Brightness	Standard Mode	80%	
				Battery Mode	20%	
		Firmware Update Info	activated			
		Sound	Volume	10%		
			Key Tone	-		
		Factory Settings		-----		
		Communication	Communication	Device name	Eurostar 60 WiCo	
				Remote Control	...	
				WiCo	activated	
Other	-					
Bluetooth	activated					
Information	Information			Stirring	Max Speed	yes
					Max Torque	yes
		Interval Run	yes			
		Interval Stop	yes			
		Safety	Time Out	yes		
Software	Software	Safe Speed	yes			
		Version	yes			
		Station Version	yes			
Firmware Update Info	Operating Mode	yes				
Firmware Update Info	Firmware Update Info	yes				

Note: ❶ Only for **EUROSTAR 100 control!**
 ❷ Only for **EUROSTAR 400 control!**

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IKA®-Werke GmbH & Co.KG