## **HG-15A**

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# Contents

<b>1 Introduction .....</b>	<b>8</b>
<b>2 Safety Instructions .....</b>	<b>8</b>
<b>2.1 Description of symbols.....</b>	<b>8</b>
<b>2.2 Instructions .....</b>	<b>9</b>
<b>3 Product description .....</b>	<b>10</b>
<b>4 Packing contents.....</b>	<b>10</b>
<b>5 Positioning the device and commissioning.....</b>	<b>10</b>
<b>6 Job description .....</b>	<b>11</b>
<b>7 Maintenance and cleaning.....</b>	<b>11</b>
<b>8 Technical data.....</b>	<b>11</b>
<b>9 Certificate.....</b>	<b>12</b>

# 1 Introduction

Thank you for purchasing our **HG-15A Homogenizer**.

You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**Risk of injury**

Avoid injuries of your hands by following these instructions.

## 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty and may cause fatal injuries.

- Do not use this device in area with potentially explosive atmosphere.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.
- Dry run may destroy the dispersing tool.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.

- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not put your hands into the device before the motor has stopped.
- Turn the device off before change dispersing tools.
- Demount dispersing tools before cleaning.

### 3 Product description

The **Wisdom Homogenizer HG-15A** is designed for nearly all media and dispersing processes in chemistry, biology, medical and pharmaceutical science.

Key features are as follows:

- Simplest operating.
- Fast commissioning.
- Long life by simply and solid construction.
- Over temperature and over current protection.
- Up to 27.000 rotations per minute.
- The dispersing tools are made of stainless steel and PTFE and are suited for all sterilization methods.

### 4 Packing contents

Quantity	Denomination		
	DH.WHG02016	DH.WHG02017	DH.WHG02015
1	<b>Homogenizer HG-15A</b>		
1	Allen key		
1	body holder		
1	operating manual		
1	ST110 standard plate stand		
1	CL200 clamp		
1	HT1018 dispersing tool	HT1025 dispersing tool	

### 5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Remove perhaps soiling and package material.
5. Attach the clamp to the tripod rod.
6. Mount the body holder at the clamp.

7. Stick the device at the body holder.
8. Tighten the bolts with the included Allen key carefully.
9. Stick the dispersing tool into the device and tighten the screw.
10. Check that each connection fits close.
11. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
12. Make sure that you follow all safety instructions.
13. Connect the device with the power outlet.
14. Switch the device on.
15. Set your values and start the process (look at point 6).

**Never run the device without any load and dip it into the medium which should be homogenized before powering on the device. Overspeed and resulting vibrations can damage the device.**

**Don't push the homogenizer constantly against the vessel, if you are homogenizing inside a glass vessel. High frequent vibrations could damage your glass vessel.**

## 6 Job description

Switch the device on and turn the knob to set the speed.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Technical data

		HG-15A
<b>Dimension (mm) (wxdxh)</b>		60x70x215
<b>Speed (U/min)</b>		2000 - 27.000
<b>Dispersing capacity</b>		up to 2,5 liters
<b>Motor rating input /output (W)</b>		300 / 160
<b>Safety device</b>		overheat protection
<b>Material</b>	<b>Housing</b>	powder coated steel
	<b>tools</b>	stainless steel and PTFE (bearing)
<b>Weight (kg)</b>		2,5
<b>Usable tools</b>		<b>HT1008, HT1010, HT1018, HT1025</b>
<b>Power supply</b>		AC 230 V, 1N~, 50/60 Hz

## 9 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

<b>Article</b>	<b>HG-15A Homogenizer</b>
<b>Serialno.</b>	
<b>Date</b>	

**witeg Labortechnik GmbH**

Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)